# Operating Instructions 使用说明书

8-Channel Adjustable Volume Pipette

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MP10 / MP50 / MP300

# Contents

1.Product Introduction	2
1.1 Product features	2
1.2 Box contents	2
1.3 Product structure	3
1.4 Specifications	3
2.Operation	4
2.1 Setting volume	4
2.2 Attaching and ejecting tips	4
3. Pipetting	5
3.1 Notes	5
3.2 Forward pipetting	5
3.3 Reverse pipetting	7
4.Performance Test and Calibration	8
4.1 Performance test	8
4.2 Calibration	8
5.Maintenance	10
5.1 Cleaning	10
5.2 Maintenance	10
6.Pipette Tips Compability	14
7.Spare Parts List	15
8.Troubleshooting	16
9.Warranty	17

# 1. Product Introduction

## 1.1 Product features


- Ergonomic design allows comfortable grip and fatigue-free pipetting
- Wide volume range: from 0.5~300ul
- 4-digit volume display provides accurate volume setting and visibility
- Easy to disassemble the lower part for replacement or cleaning
- Tip contact parts are chemical-resistant
- Easy to calibrate and maintain with tools supplied.
- Different volume range is identified by different color top
- Each pipette is calibrated in accordance with ISO8665
- The lower part can be autoclaved

## 1.2 Box contents

Item	Qty.	Item	Qty.
Pipette	1	Tip	8
Operating instruction	1	Certificate	1
Calibration tool	1	"ADJ" sticker	1
Grease	1		

### 1.3 Product structure

- ① Control button (ASA/PC)
- ② Volume setting ring (ASA/PC)
- ③ Ejector button (ASA)
- ④ Display window (PC)
- ⑤ Handle (ASA)
- ⑥ Detach button (PEI)
- ⑦ 8-Channel lower part (PEI)
- ⑧ Ejector sleeve (PEI)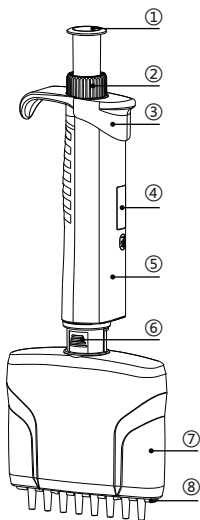


### 1.4 Specifications

Model	Volume Range ( $\mu$ l)	Testing Volume ( $\mu$ l)	Maximum permissible systematic error		Maximum permissible random error	
MP10	0.5-10	10	2.00%	0.200ul	1.00%	0.100ul
		5	4.00%	0.200ul	2.00%	0.100ul
		1	8.00%	0.080ul	5.00%	0.050ul
MP50	5-50	50	1.00%	0.500ul	0.40%	0.200ul
		25	1.50%	0.375ul	0.80%	0.200ul
		5	3.00%	0.150ul	2.00%	0.100ul
MP300	30-300	300	0.60%	1.800ul	0.30%	0.900 $\mu$ l
		150	1.00%	1.500ul	0.50%	0.750ul
		50	3.00%	1.500ul	1.00%	0.500ul

## 2.Operation



### 2.1 Setting volume

- Turn the volume setting ring shown as above to set the volume.
- The height of the control button changes as the volume is adjusted. The higher the button is, the bigger the volume will be. **DO NOT** turn the volume setting ring excessively to set the volume outside of range, which would jam the mechanism and damage the pipette.
- The side viewing window displays the set volume.
- The numbers on the volume display are to be read from the top to the bottom. The decimal places start after the underscored digit.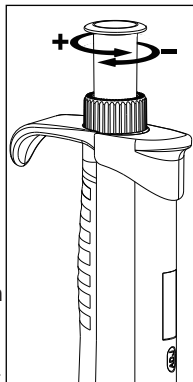

### 2.2 Attaching and ejecting tips

- The pipette can only be used after a pipette tip has been attached.
- Make sure the tip cone is clean before attaching the tip.
- Make sure it is a tight fit between the tip and the tip cone for accurate transferring of liquids.
- When attaching a tip by hand, it must be handled in a way that prevents both the contamination and hand-warming of the pipette tip.
- Press the ejector button firmly downwards to eject the tips.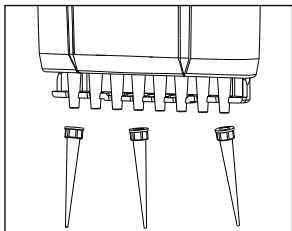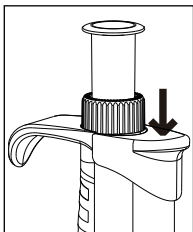

## 3. Pipetting

### 3.1 Notes


- To ensure maximum precision and accuracy, we recommend initially wetting each new tip by aspirating and dispensing the liquid three to five times.
- To aspirate liquid, immerse the pipette tip vertically in the liquid at a specific immersion depth (refer table below) and maintain the immersion depth to ensure no air is aspirated. Carefully wipe the pipette tip against the tube inner wall to ensure that no outer wetting remains on the tip.
- To dispense liquid, place the tip on the tube inner wall at an angle. Hold down the control button to dispense the liquid and wipe the tip on the tube inner wall.

### 3.2 Forward pipetting


- Press down the control button to the first stop.
- Vertically immerse the pipette tip into the liquid at a specific immersion depth (refer table below)

Sample volume	Immersion depth	Waiting time
0.5-100ul	2-3mm	1s
100ul-300ul	2-4mm	1s
> 1000ul	3-6mm	3s


- Allow the control button to slide back slowly. Maintain the immersion depth to ensure no air is aspirated.
- Wait before removing the pipette tip from the liquid. The waiting time is based on how large volume is aspirated.
- To dispense the liquid, press down the control button slowly until the first stop.
- To empty the tip, press the control button until the second stop.
- Allow the control button slide back to the starting point slowly.
- To eject the tip, press the ejector button.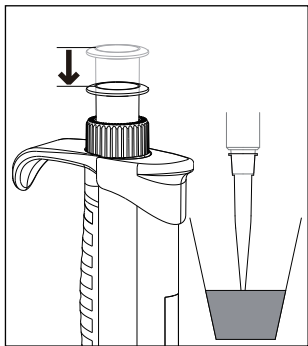


1. Press down the control button to the first stop.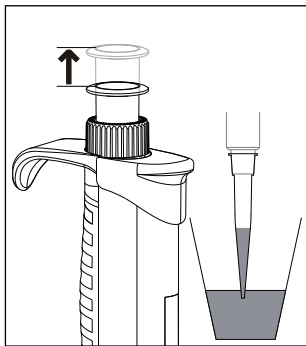

2. Let the control button slide back slowly.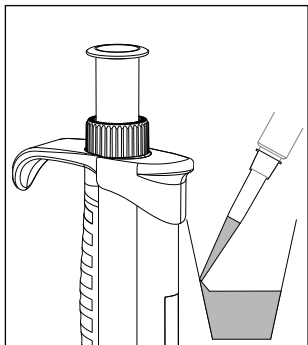

3. Place the pipette tip against the tube inner wall, at an angle of around 45°. Press the control button slowly to the first stop position and wait until the flow of liquid stops.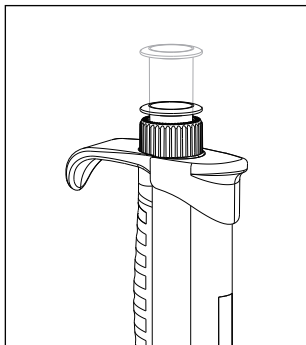

4. To empty the tip, press down the control button to the second stop. Hold down the control button and wipe the tip on the tube inner wall.

### 3.3 Reverse pipetting

Reverse pipetting is used to dispense bubbling liquid, highly viscous liquid and can also be used for very small volume. Before pipetting, please wet the tip by aspirating and dispensing the liquid.

- Press down the control button to the second stop.
- Vertically immerse the pipette tip approx. 2-3mm into the liquid.
- Allow the control button to slide back slowly. Maintain the immersion depth to ensure no air is aspirated.
- To dispense the liquid, press the control button slowly until the first stop.
- The remaining liquid together with tips should be discarded.

## 4. Performance Test and Calibration

### 4.1 Performance test

Performance test should be performed in a draught-free room at 15-30°C (constant to  $\pm 0.5^\circ\text{C}$ ) and the relative humidity above 50%. Keep the pipette, tips and test water in the test room for at least 2 hours prior to the test. Please perform the tests with the distilled water or deionized water (grade 3, ISO3696) and the analytical balance (accuracy up to 0.01mg).

- Attach the tip to the pipette properly.
- Wet the tip by aspirating and dispensing the set volume 3 to 5 times.
- Press down the control button to the first stop
- Vertically immerse the pipette tip approx. 4mm into the liquid.
- Allow the control button to slide back slowly. Maintain the immerse depth to ensure no air is aspirated.
- Dispense the liquid into the weighing container and place the tip on the container inner wall at an angle. Hold down the control button and wipe the tip on the container inner wall.
- Read the weight in mg.
- Repeat the same tests for 10 times and record each reading.
- Convert each weight reading to volume by multiplying the error correction Z ( $Z=1.0033$ ,  $22^\circ\text{C}$  and  $101.3\text{ Kpa}$ ).

### 4.2 Calibration



- Remove the "ADJ" sticker.
- Insert the calibration tool into the hole where the "ADJ" sticker is

located. Please make sure the wedge portion is facing downwards.

- Turn the volume setting ring to adjust the volume. To increase the volume, adjust in the counterclockwise direction and to decrease the volume adjust in clockwise direction.
- Rotate the volume setting ring counterclockwise to increase the displayed volume (if it is smaller than the actual volume) or rotate the volume setting ring clockwise to decrease the displayed volume (if it is larger than the actual volume) until the displayed volume is same with the actual volume.
- Attach the "ADJ" sticker after calibration.
- Perform tests as mentioned above to verify the calibration.

**Note:** Each pipette is calibrated at our facility before shipment.

Recalibration is to eliminate the errors caused after long-term use.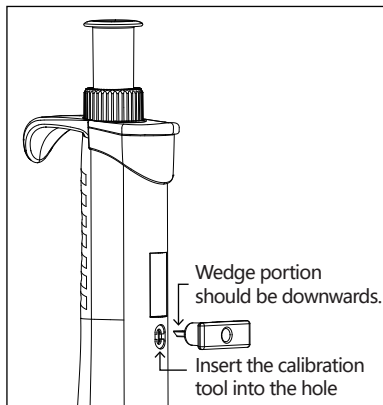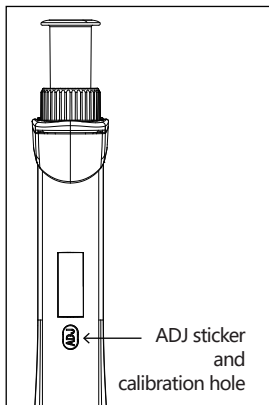

## 5. Maintenance

Aggressive substances may damage the pipette, pipette tip and accessories. Faulty dispensing results are sometimes due to lack of maintenance.

- Check the resistance to chemicals before using organic solvents or aggressive chemicals.
- Only use liquids whose vapors do not corrode the used materials.


### 5.1 Cleaning

The lower parts are wear parts. Clean them after contamination or use of aggressive chemicals. If the lower parts are worn or damaged, replace these parts. Unsuitable cleaning agents can damage the device.


- Moisten a cloth with a cleaning agent / alcohol.
- Remove external contamination.
- Moisten a new cloth with water.
- Wipe down the housing.

### 5.2 Maintenance


- Press both sides of the ejector sleeve to detach the lower part.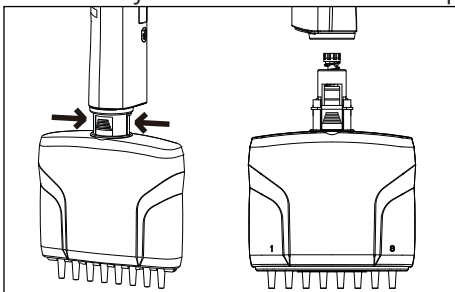


- Slide the cover plate downwards to open (The side with number "1" & "8" should face the operator).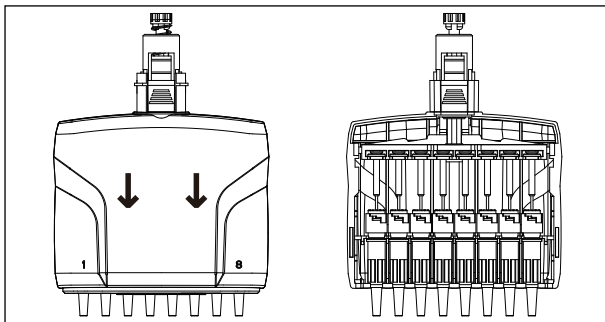


- Press down and rotate to take out the spring mount and the spring.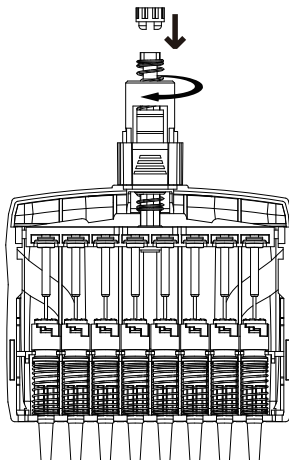

- Push down the slide block towards the piston cylinders. After this, carefully release the piston cylinder. Insert a tool (like screwdriver) below the top portion of the piston cylinder and move it upward to release it from its position. ( To reinsert the piston cylinder, the bottom portion should be secured in the slot first, and then push down the top into position.)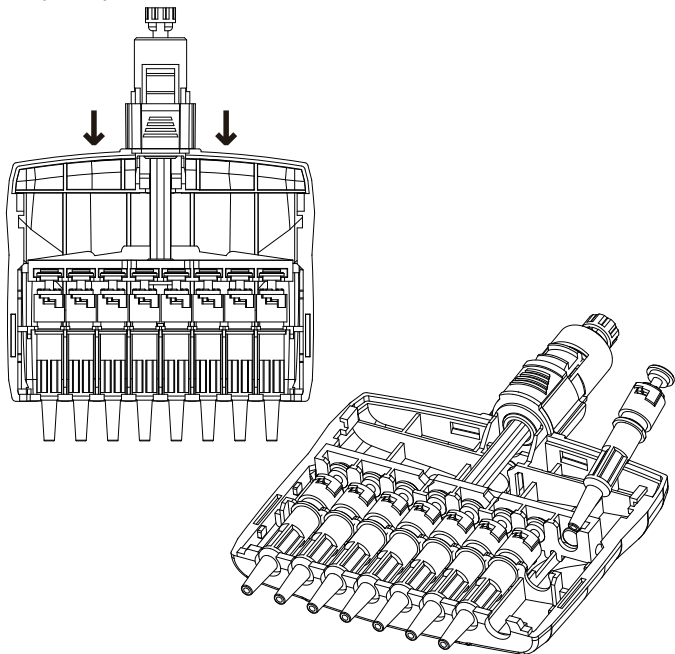

- To disassemble the piston cylinder, press down and rotate the piston mount to separate the piston spring, O-ring and piston guide.

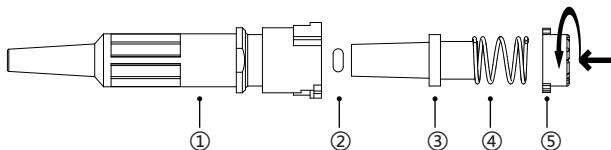- ① Piston cylinder
- ② O-ring
- ③ Piston guide
- ④ Piston spring
- ⑤ Piston mount

- Wipe the piston cylinder, piston pin, spring, fixed seat, O-ring and ejector sleeve. After cleaning, brush moderate amount of silicone grease onto the piston pin, fixed seat and O-ring.

**Note:** excessive silicone grease may interfere with smooth movement of the piston.

- After re-assembly, push the control button several times to make sure the silicone grease is uniform. (The pipette is without any liquid when doing this operation.)

## 6. Pipette Tips Compatibility

Brand	Volume	MP10	MP50	MP300
Eppendorf	10uL			
	200uL			
	300uL			
Axyen	10uL(universal)			
	10uL(long)			
	200uL			
	300uL			
QSP (Thermo scientific)	10uL			
	200uL			
Gilson	10uL			
	200uL			
Kirgen	KG1011(0.1-10uL)			
	KG1111(0.5-10uL)			
	KG1212(1-200uL)			
	KG1213(1-300uL)			
Cellpro	10uL			
	200uL			
	300uL			
Pakgent	UFRT-10R			
	UFRT-10R-L			
	UFRT-10ER			
	UFRT-10ER-L			
	UFRT-200R			
	UFRT-200R-L			
	UFRT-200ER			
	UFRT-200ER-L			

## 7. Spare Parts List

	Part No#			
Volume Range	Piston cylinder	Piston rod	Tip O-ring	Piston spring
0.5-10 $\mu$ l	3.1.305235000	2.3.205235100	/	3.2.405222004
5-50 $\mu$ l	3.1.305227000	2.3.205227100	3.4.605222001	3.2.405222004
30-300 $\mu$ l	3.1.305222009	2.3.205222100	3.4.605222001	3.2.405222004

## 8. Troubleshooting

Problem	Cause	Solution
Residual liquid and/or liquid is dripping from the tip and /or the dispensed volume is too small.	<ul style="list-style-type: none"> <li>• Incompatible tip.</li> <li>• The tip is not wet uniformly.</li> <li>• Tip incorrectly attached.</li> <li>• Foreign particles between the tip and the piston cylinder.</li> <li>• Insufficient grease on the piston and O-ring; O-ring positioned incorrectly, or the O-ring is damaged.</li> <li>• Malperation; The calibration is changed, or the pipette is damaged.</li> </ul>	<ul style="list-style-type: none"> <li>• Use compatible tip.</li> <li>• Attach new tip.</li> <li>• Attach the tip firmly.</li> <li>• Clean the piston cylinder and attach new tip.</li> <li>• Clean and grease the O-ring and piston. Reinstall or replace the O-ring.</li> <li>• Recalibrate and repair the pipette</li> </ul>
The control button jams and does not move smoothly.	The piston is dirty. The pipette is blocked.	Clean the piston cylinder and ejector sleeve.
The pipette is blocked, or the dispensed volume is too small.	Insufficient silicone grease on O-ring. Foreign particles on the piston pin and O-ring.	Clean and lubricate the O-ring and piston. Clean the end of piston cylinder.
The ejector sleeve is blocked or does not move smoothly.	The spring jams.	Grease the spring.

## **9.Warranty**

Each pipette is calibrated and tested before shipping and it's ready for use.

The pipette is covered by warranty for one year against defects in materials and workmanship. If your pipette failed to function, please contact your local representative.

Any warranty will, however, be deemed as void if fault is found to have been caused by maltreatment, misuse, unauthorized maintenance of service or negligence of regular maintenance and service, accidental damage, incorrect storage or use of the products for operations outside their specifications, contrary to the instructions given in this manual or with other than the manufacturer's original tips or compatible tips recommended by the manufacturer.

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